

Installing ACD spectrus for basic data processing

The University has a site license for ACD Spectrus Processor (running under Windows), which is a piece of software that can be used for processing NMR data. To install the software the following steps should be taken.

1. In **File Explorer** or **My Computer** navigate to
\\ss7a.ds.man.ac.uk\vol5\vol3\users\snmrdata\ACD_spectrus_processor_download or
\\nasr.man.ac.uk\epsr\$\$\snapped\replicated\NMR_ACD_spectrus_processor_download
2. Double-click **ACDLabs2015p65_SpectrusProcessor_LSIInstall.exe**.
3. If you get a 'security pop-up' click **Yes**.
4. Click **Next**. Then read and accept the terms in the License Agreement, and click **Next**.
5. Enter **10.99.106.130** as server address and **7199** as the port number.
6. Deselect any components of the ACD software that you do not wish to install, then click **Next**.
7. Choose a folder then click **Next**, or just click **Next** to proceed with the default folder.
8. Deselect *Install ChemBasic Goodies* unless you want it installed. Click **Next**.
9. Click **Install**. Wait for the software to be installed. It will also update to the latest version.
10. Click **Finish >>**.
11. Click **OK** to restart the computer and complete installation.
12. If you are asked for a web update key, enter: **Q419QNU7EWFHWMWQ8K0P**.
13. To get you started read the quick start guides in the
\\ss7a.ds.man.ac.uk\vol5\vol3\users\snmrdata\ACD_spectrus_processor_download
directory.

The use of ACD spectrus processor requires connection to the University of Manchester network. If you wish to use it off campus please connect via VPN. Instructions for connecting via VPN can be found on the IT services web pages: <http://www.itservices.manchester.ac.uk/our-services/my-it/vpn>.